*Fs flame*¹⁰⁰ *Safety Bunsen*

Sicherheitsbunsenbrenner
Safety Bunsen Burner

Windschutz / Windshield
Edelstahl / stainless steel
Art.-Nr./Art.-No. 2.000.600


Spritz- und Wachsschutz
Edelstahlplatte und Glaszylinder
Spray and Wax protector
stainless steel base and glass cylinder
Art.-Nr./Art.-No. 2.000.200
ohne Glaszylinder / w/o glass cylinder
Art.-Nr./Art.-No. 2.000.201

Externer DoubleClick IR-Sensor
zum berührungslosen Flammenstart
External DoubleClick IR-Sensor
for touch-free flame activation
Art.-Nr./Art.-No. 6.000.406

(1) Fußpedal / Foot pedal
Edelstahl / stainless steel
Art.-Nr./Art.-No. 6.000.402
(2) Fußpedal Mini / Foot pedal Mini
Kunststoff / plastic
Art.-Nr./Art.-No. 6.000.403
(3) Tischtaster / Benchtop switch
Edelstahl / stainless steel
Art.-Nr./Art.-No. 6.000.405
(4) RF-Stick Funkset / RF-Stick radio set
kabelloses Edelstahl-Funkfußpedal
stainless steel radio foot pedal
(available in EU countries only)
Art.-Nr./Art.-No. 2.000.404-RF

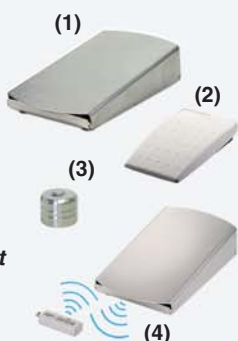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* optional

Flame¹⁰⁰ – sets new standards in the Bunsen burner class!

The safe alternative for all traditional Bunsen or alcohol burners!

Multi-Use

The Flame¹⁰⁰ is ideally suited for all flame related applications in the laboratory. The precision flame allows graduated heating of dental tools as well as safe sterilization of microbiological instruments. Flame size and intensity can be adjusted infinitely.

Easy operation

The safety Bunsen burner Flame¹⁰⁰ activates immediately with the push of a button. No match or a pilot flame are required. Optionally, the Flame¹⁰⁰ can also be operated by a foot pedal* or an external infrared motion sensor*. The Flame¹⁰⁰ is suitable for stationary natural gas and propane / butane gas supplies as well as gas cartridges or gas cylinders. Numerous adapter systems are optionally available.

Reliable safety

The proven Safety Control System (SCS) is also incorporated in the Flame¹⁰⁰. All potential hazards are constantly monitored and, if necessary, protective measures – such as the shutting off the gas supply – are activated. In addition to the ignition and flame control function and the overheating protection the safety package also features the continuous burner head control BHC. BHC unfailingly identifies burner head clogging by liquids or solid substances. Compared with a conventional Bunsen burner, the Flame¹⁰⁰ with the safety standard „SCS“ eliminates the risk of a gas leak or an explosion.

Efficient

The Flame¹⁰⁰ is economical to operate and attractively priced for any laboratory, thus making savings for both your budget and the environment. Compared with traditional Bunsen burners, the gas consumption is reduced to a minimum, because the flame burns only when it is really needed!

Compact and rugged

The housing and burner tube of the Flame¹⁰⁰ are made of stainless steel. The removable burner tube makes cleaning easy and the small foot print is an advantage for any lab bench.

Flame¹⁰⁰ - for a safe flame in the laboratory!

Neigungsverstellung / Tilt adjustment
Edelstahl / stainless steel
Art.-Nr./Art.-No. 2.000.300

Tray / Tray
Edelstahl, 5 Ablagemöglichkeiten
stainless steel, 5 places
Art.-Nr./Art.-No. 6.000.320

Adapter 360 / Adapter 360
für Kartuschen CV 360
for gas cartridges CV 360
Art.-Nr./Art.-No. 2.000.701

Programs

Button function	✓
Pedal Standard (foot pedal optional)	✓
Safety Control System SCS®	✓
Burner Head Control BHC	✓
Gas inlet filter	✓
Removable & decomposable burner head	✓
Housing, burner head, gas connection & controls made of stainless steel	✓
DVGW approved	✓

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Art.-No. 2.100.000

with button function

Safety Control System (SCS) with Burner Head Control (BHC)

2 standard programs for button (function knob) and foot pedal (foot pedal optional)

Removable burner tube

Holding device for 2 inoculation loop holders

Nozzles for natural gas, propane / butane gas

Wrench 17 mm for gas connection

Tubing connector with swivel nut

Screwdriver for electrode holder

Switching power supply (global)

Instruction manual and 2-year warranty

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Start-Stop with timer, 60 min

Foot pedal:

Standard (flame during pressed foot pedal, foot pedal optional)

Safety features

Safety Control System (SCS)

with gas safety cut off:

ignition and flame control
temperature monitor,
burner head clogging monitor (BHC)
automatic unit switch off, 4 h

Gas supply and consumption

Gas connection:

1/4" left with gas filter

Gas types:

II2ELL3B/P: natural gas E/LL, 18 - 25 mbar
liquid gas, 20 - 50 mbar

Connected load:

48 g/h liquid gas, 53 l/h natural gas (LL)

Continuous cartridge operation:
(approx.)

CV 360 - 65 min, CG 1750 - 210 min
C 206 - 230 min, CP 250 - 305 min
CV 470 - 550 min, CV 300 - 290 min

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WLD-TEC GmbH